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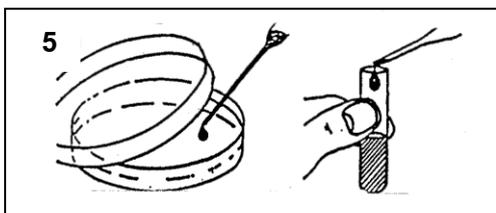
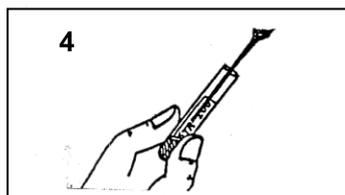
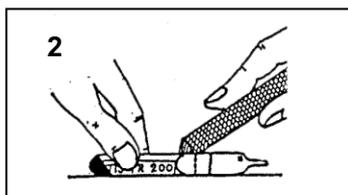
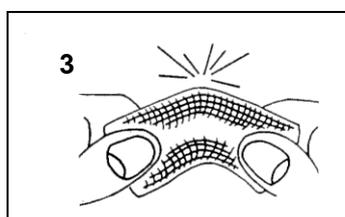
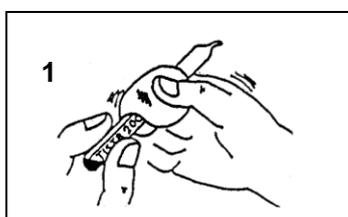
Distribution Form

SENT TO :
.....
.....

Table with 6 columns: TISTR No., Name of Microorganism, Preservation Date, Best before (D/M/Y), Medium, Conditions. The table contains 10 empty rows for data entry.

- Note : 1. Ampoules have not opened soon after receipt should be stored in the dark, preferably in a cool place.
2. Third party distribution is prohibit.
3. TISTR Culture Collection kindly requests the customers to forward the strain information/ publications obtained from further investigation of the strains.

DEBIT NOTE No. :
PACKED BY :
DATE :

Thailand Institute of Scientific and Technological Research (TISTR)

35 Moo 3 Technopolis, Klong 5, Klong Luang,

Pathum Thani 12120, Thailand

Tel: (662) 577 9034, (662) 577 9041 Fax: (662) 577 9031

E-mail: mircen@tistr.or.th

Revival of Lyophilized or Freeze-Dried Cultures

1. After sterilizing ampoule with 70% alcohol-dampened gauze, make a file cut on the ampoule at the mid-point of the cotton wool plug.
2. Wipe the ampoule with 70% alcohol-dampened gauze.
3. Carefully break the ampoule at the scored area, break it gently.
4. Aseptically add to the freeze-dried culture 0.3 to 0.4 ml of an appropriate liquid medium with a Pasteur pipette; mix well.
5. Transfer a one drop of the suspension and streak on to the surface of an appropriate agar plate and rest of the suspension to a test tube containing 5 ml of the same liquid medium. [For fungi will be transfer three drops (make three points) of the suspension on to the surface of an appropriate agar plate just drop not streak]
6. Incubate the culture (Both agar plate and liquid medium) at the appropriate temperature and allow sufficient time for growth to occur.